

Adjustable Film Applicator

BEVS 1806

User Manual



Version 201412

This manual shall be read carefully before starting. Directions included in this operation manual shall be strictly followed.

4. Calibration

Before use check the zero settings of the micrometer heads as follows:

1. Place applicator on a polished plate glass of suitable size. Set micrometer heads to zero.
2. When the applicator is placed between the observer and a bright light the blade should touch the glass from end to end and not lift the end section. The zero point should check within ± 1 , division of the micrometer - this is set at manufacture.

If applicator does not zero and resettings are required proceed as follows:

Set the micrometer heads to zero, loosen the micrometer clamp set screws, press the knife blade down onto the glass plate, press down the micrometer heads until their spindles are contacting the blade and then tighten the micrometer clamp screws.

5. Maintenance

1. There will have the film error as the knife blade appears crack and transformation by the falling and crash. So take good care of the equipment when it is in process of use.
2. Screw is not allowed to tear down as the calibration was already finished before delivery. Or it will have the error.

6. Order information

BEVS 1806/50	Adjustable Applicator	50mm Width
BEVS 1806/100	Adjustable Applicator	100mm Width
BEVS 1806/150	Adjustable Applicator	150mm Width
BEVS 1806/200	Adjustable Applicator	200mm Width
BEVS 1806/250	Adjustable Applicator	250mm Width
BEVS 1806/300	Adjustable Applicator	300mm Width